

Double staining LYVE-1/CD31 on humanen cryo-sections

- dry cryo sections for 1h at RT
- fix sections for 10 min at -20°C in MeOH
- wash 3x3 min in PBS at RT
- **block 20 min in 10% goat serum in PBS**
- **1. AB over night at 4°C in 10% goat serum in PBS**
[anti h-LYVE-1 10µg/ml and anti h-CD31 1:50 (DAKO Klon JC70A Code No. M0823)]
- wash 3x3 min in PBS at RT
- **2. AB 30 min at RT in PBS**
[goat anti rabbit IgG (H+L) CY3 1:500 (dianova 111-165-144) and goat anti mouse biotin 1:500 (DAKO Code No. E0433)]
- wash 3x3 min in PBS at RT
- **staining of nucleus 10 min at RT in Hoechst Dye 1:5000 in PBS**
- wash 3x3 min in PBS at RT
- **cover with DAKO Fluorescent Mountingmedium**